



Sterile ready to use solution based on peracetic acid for disinfecting surfaces in cleanrooms.

perform® sterile PAA

Our Plus

- fast, effective and sporicidal
- sterile (filled aseptically in cleanrooms A/B)
- microbial filtered (0,2 µm)
- ready to use in a handy, sterile 500 ml spray system
- colour and perfume free
- efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing Practice: Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing Practice; U.S. Department of Health and Human Services, Food and Drug Administration, September 2004.

Instructions for use

Apply perform® sterile PAA undiluted to pre-cleaned objects and surfaces and allow to take effect, ensuring that all surfaces are moistened completely.

Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal EN1276, EN13697	ready-to-use	5 min.
fungicidal EN1650, EN13697	ready-to-use	5 min.

Efficacy	Concentration	Contact time
virucidal against enveloped viruses in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline	ready-to-use	1 min.
sporicidal EN13697, EN13704	ready-to-use	5 min.



Product data

Composition:
100 g solution contains the following active substances: 0.09 g peracetic acid.

Chemical-physical data

Color	colourless
Density	1,01 g/cm ³ / 20 °C
Flash point	> 105 °C
Form	liquid
pH	ca. 3,5 / 10 g/l / 20 °C / in water
Viscosity, dynamic	1,05 mPa*s / 20 °C

Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: www.schuelke.com. For individual questions:

Customer Sales Service
Phone: +49 40 52100-666
E-Mail: info@schuelke.com

Manufacturer:
BIOXAL SA
Route de Varennes
BP 30072
71103 Chalon sur Saône
France

Special advice

Use disinfectants safely. Always read the label and product information before use.

Professional use only. Read safety data sheet before use.
Good material compatibility with stainless steel, PTFE, PVDF, PEHD, PP, Viton, glass, etc.
Contact with non-ferrous metals (brass, copper) should be avoided. Wearing of protective gloves (e.g. made of Nitrile rubber) is recommended. Carry and store product container in an upright position.

Information for order

Item	Delivery form	Item no.
perform [®] sterile PAA 500 ml bottle	10/Carton	on request
perform sterile PAA 300 ml FL	10/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

Application aids

Application aids	Item no.
perform [®] spray bottle holder	693540
perform [®] spray bottle holder - plate	693541

Related Products

- perform[®] sterile dry wipes
- perform[®] sterile mix dry wipes
- perform[®] sterile premium dry wipes

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining our high quality standards.



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com